

## Micro Learning Video: CCC Confirmation of Interpretation Factors

In this video, you will learn how to review and verify the interpretation factors of a specimen.

1. Click the Accession # in the Specimens list view to open a specimen record.
2. Click the Pregnancy Information tab.
3. Next, validate the Patient Interpretation Factors.  
Click the checkbox next to each field to validate it.
4. Next, validate the Case Interpretation Factors.  
Click the checkbox next to each field to validate it.
5. Next, validate the Specimen Interpretation Factors.  
Click the checkbox next to each field to validate it
6. Click Save.
7. Alternatively, click Save & Close.